



BioPAT® Xgas

Online Off-Gas Analysis for Monitoring and Controlling Bioprocesses

Benefits

- One-sensor determination of O₂ and CO₂
- Integration into BIOSTAT® and BioPAT® MFCS
- Monitoring of metabolic processes
- Process and yield optimizing tool
- Compact design for flexible use
- Automatic moisture and pressure compensation



Product Information

The BioPAT® Xgas, can monitor exhaust gas from bioprocesses in real time. It is a dual oxygen-carbon dioxide sensor featuring automatic compensation of humidity and pressure. The oxygen concentration can be measured in a range of 1 – 50% and the carbon dioxide concentration in a range of 0 – 10%.

O₂ Sensor

The zirconium-dioxide sensor measures the oxygen concentration over a range of 1 to 50% with an accuracy of under 0.2% full scale over the whole measurement range with a $\pm 3\%$ reading on the value. The measurement principle requires the sensor temperature to be 580°C (1,076°F) and operates to specification for up to 15,000 hours (approx. 20 months) with less than 2% drift. The O₂ sensor requires a minimum of 0.1% oxygen during operation to avoid damage to the electrode cell. Therefore, sensor exposure to non-oxygen gas mixtures must be avoided.

CO₂ Sensor

The dual-wavelength infrared sensor detects the carbon dioxide concentration over a range of 0 to 10% with an accuracy of under 0.2% FS with over the whole measurement range with a $\pm 3\%$ reading on the value. The robust sensor operates for up to 20,000 working hours (approx. 25 months) with less than 2% drift. As a nonstandard, the sensor can be recalibrated to measure higher concentrations of CO₂ at a lower accuracy.

Dimensions and Exhaust Connection

The BioPAT® Xgas sensor has a compact footprint and a stainless steel housing (IP65) that connects to standard gas flow path adapters. This allows the sensor to be used at different maximum gas flow rates for each size of bioreactor. Sartorius Stedim will provide the appropriately sized flow path (1/4" to 1 1/4") for each BIOSTAT® application and arrange for mechanical and software installation. The system is mountable and can be integrated into the Sartorius Stedim bioreactor vessel exhaust line with either stainless steel piping or flexible tubing. This means a BioPAT® Xgas is dedicated to a single cultivation vessel with an adjustable measurement frequency down to every 10 seconds. It is recommended to mount the BioPAT® Xgas at a downward angle with respect to the inlet gas flow direction in order to minimise water collection in the measurement flow path.

Pressure and Humidity Sensor

As pressure and moisture levels affect the detected level of the O₂ and CO₂, these levels must be measured and compensated for. Internally, the BioPAT[®] Xgas has a piezoresistive silicon pressure sensor (0.8 to 1.3 bar, 11.6 – 18.85 psi) and a capacitive polymer humidity sensor (5 to 100% RH) sensor. These sensors account for and compensate over an exhaust gas temperature range of 15 to 40°C, or 59°F to 104°F.

Data Transfer

The sensor data can be transferred to the BioPAT[®] DCU as an analog signal (4 – 20 mA) or as a voltage signal (1 – 10 V). An OPC[®] interface to our SCADA software BioPAT[®] MFCS/win allows digital data transfer. For either connection method, the data cables and digital conversion boxes are provided by Sartorius Stedim in a choice of integration packages, either for computers or for BioPAT[®] DCU.

BioPAT[®] DCU

The BioPAT[®] Xgas sensor readings are shown both on the visual display and on the batch record plot of BioPAT[®] MFCS as two separate output signals for O₂ and CO₂ exhaust percentages. A retrofit of an existing BIOSTAT[®] will require a DCU configuration update and two available external input connections.

BioPAT[®] MFCS

Regardless of whether the sensor data comes from the BioPAT[®] DCU or is transferred by OPC[®] to BioPAT[®] MFCS, it is handled in the same way. The BioPAT[®] Xgas Vis software functions as the OPC[®] server, sending data to the BioPAT[®] MFCS/win OPC[®] DA Client for up to 12 sensors. Once the data is in BioPAT[®] MFCS, it can be used in conjunction with the data from the mass flow controllers and other BIOSTAT[®] peripherals to calculate, plot and track the process metabolic and respiratory factors. These factors can be used to trigger actuators for automating gas-mixing strategies, adding feed or induction media to increase process yield and consistency.

Accessory Requirements

All power and data transfer cables are provided for the choice of connection method offered. The integrated feature for pressure and humidity compensation means no moisture traps, gas coolers or pumps are needed to operate the BioPAT[®] Xgas. For steam-in-place bioreactors (BIOSTAT[®] Cplus and D-DCU), a pair of automated check valves must be incorporated in the exhaust outlet setup in order to prevent the sensor from being exposed to high-temperature steam during the sterilization cycle. To enhance the precision of metabolic calculation, it is recommended to install mass flow controllers on the gas inlet lines. These controllers will allow optimal gassing settings and minimal fluctuation in DO control. The BioPAT[®] Xgas[®] has its own M12 8-pin power supply allowing it to maintain the measuring temperature and preventing condensate formation inside the sensor housing. A base plate is included with every Univesse[®] BioPAT[®] Xgas enabling it to be wall mounted for ergonomic improvements.

Metabolic Calculations

The BioPAT[®] Xgas sensor monitors the concentration of oxygen and carbon dioxide online, allowing real-time measurement of key parameters, such as oxygen uptake rate (OUR), carbon dioxide evolution rate (CER) and respiratory quotient (RQ). Beyond this, oxygen and carbon dioxide measurements enable mass balances to be calculated and the oxygen transfer rate (OTR) to be determined directly. For in-depth details, please see the BioPAT[®] Xgas Metabolic Calculations data sheet available from Sartorius Stedim.

Calibration and Interface

The unit's status can be read and a one-point calibration carried out using the two push buttons located directly on the BioPAT[®] Xgas. Annual service and factory recalibrations are recommended if the system is used for critical process control loops. The LEDs are color-coded, and a color-code chart is shown on the housing between the two manual input buttons. The exhaust O₂ and CO₂ sensor information can be obtained from the BioPAT[®] DCU or MFCS.

Technical specifications

Concentration ranges¹

0 – 10 vol% CO₂, 1 – 50 vol% O₂

O ₂ sensor unit	
Measured gas	O ₂
Measuring principle	Zirconium dioxide
Accuracy	< 0.2% FS ² ± 3% Rdg ² on value
Drift	< ± 2% value/year
Lifetime of sensor element	Approx. 15,000 operating hours
Temperature inside the sensor unit	580°C 1076°F

CO ₂ sensor unit	
Measured gas	CO ₂
Measuring principle, infrared	Dual wavelengths
Accuracy	< 0.2% FS ² ± 3% Rdg ² on value
Drift	< ± 2% value/year
Lifetime of optical components	Approx. 2 years
Temperature inside the sensor unit	3°C 5.4°F higher than process temperature

OPC® PC Connection box	
Number of Serial Ports	1
Connect directly using Telnet/SSH by port and IP address Connect with EasyPort menu by Telnet/SSH	
Serial Port Interface	EIA232/422/485 on DB9M, RJ45, DB25M or DB25F
Serial Port Speeds	50 bps to 230 Kbps with customizable baud rate support
Parity	Odd, Even, Mark, Space, None
Local Console	Port RS232 on Serial Port
Power Supply	120 V AC (USA), 230V AC (International) Wall Power Adaptor included
Nominal Input	Voltage 12v DC
Weight	0.23 kg (0.5 lbs)
Dimensions	91 × 64 × 24 (mm), 3.6 × 2.5 × 0.92 (in) case dimensions not including mounting tabs, 91 × 89 × 24 (mm), 3.6 × 3.5 × 0.92 (in) includes mounting tabs

General	
Temperature range	15 – 40°C 59°F – 104°F
Pressure range	0.8 – 1.3 bar 11.6 – 18.85 psi absolute pressure
Operating humidity	5 – 100% RH, integrated humidity compensation
Housing, stainless steel	IP65 protection
Dimensions and weight	W × L × H = 170 × 150 × 120 mm (6.69" × 5.91" × 4.72") ³ weight = 4 kg (8.82 lb)
Mechanical connection	4 mm – 1¼" ¹
Materials in contact with exhaust gas	Stainless steel, thermo-elastomer, sapphire, PTFE, polymer H.L., nitrile
Filters	PTFE, 0.22 µm; PTFE, 5 µm
Power supply	24V 1A
Storage conditions	0 – 60°C (32 – 140°F); 5 – 75% RH (non-condensing)




Electronic connections	
Power supply	8-pin M12, male
Output connection	8-pin M12, female
Electronic output	Active output, maximum 500 ohms with a 24 V power supply RS232, RS485 Modbus, 2×4 – 20 mA, USB, Modbus OPC® Server
Maintenance	One-point calibration with ambient air (0.04 vol % CO ₂ , 20.97 vol % O ₂) once a month (other conditions possible); optional factory calibration once a year
CE/FCC/ICES	EN 61326-1:2006 FCC 15:2009 Subpart 107/109, ICES-001:2006
Remarks	High concentration of NH ₃ or O ₃ could minimize the lifetime of the O ₂ sensor element

¹ Expanded ranges on request

² FS = full scale, Rdg = measured accuracy

³ Dependent on flow adapter

Technical Data

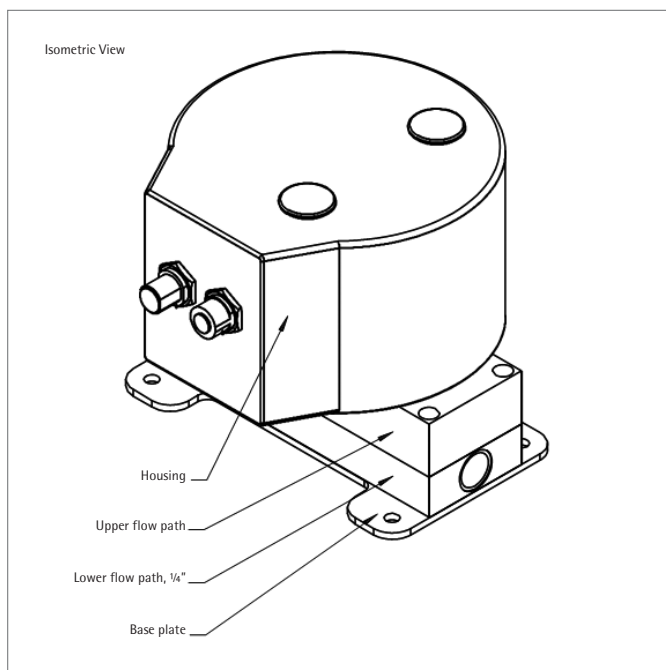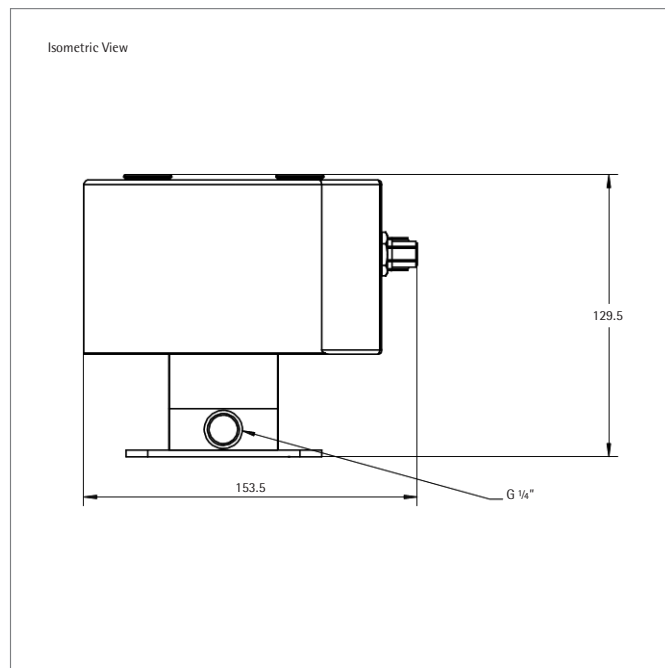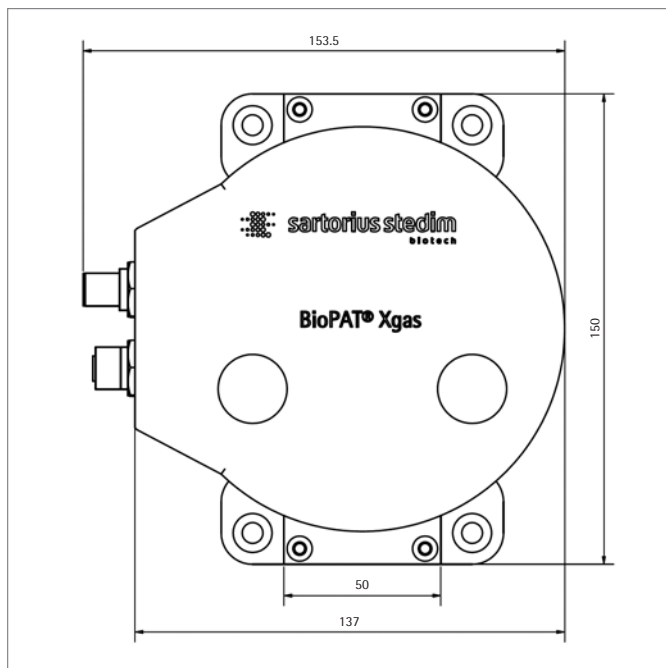

All dimensions given in millimeters

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany
Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

USA Toll-Free +1.800.368.7178
UK +44.1372.737159
France +33.442.845600
Italy +39.055.63.40.41
Spain +34.90.2110935
Russian Federation +7.812.327.5.327
Japan +81.3.4331.4300
China +86.21.68782300

Specifications subject to change
without notice. Printed and copyrighted
by Sartorius Stedim Biotech GmbH. | W
Publication No.: SLL2007-e141104
Order No.: 85032-543-77
Ver. 11 | 2014